

# Encapsulation Unit – Var J1



**Nisco Engineering AG**  
Wehntalerstrasse 562  
CH-8046 Zurich, Switzerland


Tel: +41 44 380 06 30  
Fax: +41 44 380 06 31  
e-mail: [mailbox@nisco.ch](mailto:mailbox@nisco.ch)  
<http://www.nisco.ch>

In most applications involving immobilisation of living cells or other biological materials the bead size is needed to be small (<1 mm) and carefully controlled. The reason for this is mostly because of diffusion limitation of nutrients within the hydrogel beads. An easy way for production of small alginate beads in a controllable manner is the use of a coaxial bead generator.

The basic principle of the instrument is the use of a coaxial air stream to pull droplets from a needle tip into the gelling bath. The Nisco J1 is designed for production of smaller quantities of spherical alginate beads ranging in size down to around 500  $\mu$ m.

The bead generator with coaxial airflow is basically made of Polyetheretherketone (PEEK) and can withstand most chemicals and high temperatures, which makes the unit very suitable for any disinfecting/cleaning method, typically autoclave cleaning.

*under license of FMC Polymer a.s., Oslo, Norway*

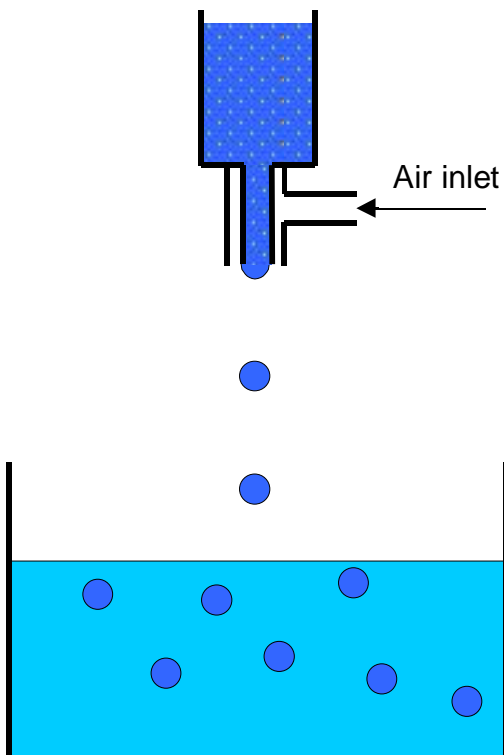

Principle of coaxial airflow bead generator



The Unit is equipped with two connections:  
. one for the hose, which feeds the alginate (or other) solution,  
. and the second connection which is meant for an air-hose with 4 mm OD.

The alginate (or other) solution may be fed into the unit with a syringe, using a syringe pump.

The magnetic stirrer is placed underneath the gelling bath to keep the beads separated during gelling.