

Heating Chamber designed for Unit – VARD2Go VAR D classic



Nisco Engineering AG
 Wehntalerstrasse 562
 CH-8046 Zurich, Switzerland

Tel: +41 44 380 06 30
 Fax: +41 44 380 06 31
 e-mail: mailbox@nisco.ch
<http://www.nisco.ch>

Laminar jet break up has some limitations concerning viscosity. Some matrices have lower viscosity with increasing temperature. Thus it is helpful for the process to work with an increased temperature. Some low melting matrices such as agarose or gelatine have to be kept in a melted condition until they have passed the nozzle. In these cases temperatures up to 60°C are sufficient. For this reason Nisco Engineering AG has developed a heating chamber in which only the syringe pump is placed. The operating display is outside the heating chamber, so that the syringes can always be kept warm.

It is important to use a syringe pump which is designed for the working temperature. We have chosen the KD Scientific pump of the 200 Series.



On the picture above you see the contained vessel of VAR A with the heating chamber.


The heating chamber is controlled by a digital PID Controller and a pT-100. The unit is delivered in the following two versions:

For 110 V

For 230 V

The chamber is designed for the following syringe pump KDS 200:

The syringe pump is controlled by a micro controller and suited to work with two parallel syringes. The specific parameters can be configured by means of display and keyboard on the pump itself. There is a TTL and a RS232 interface on the pump. The most important technical data are listed below:

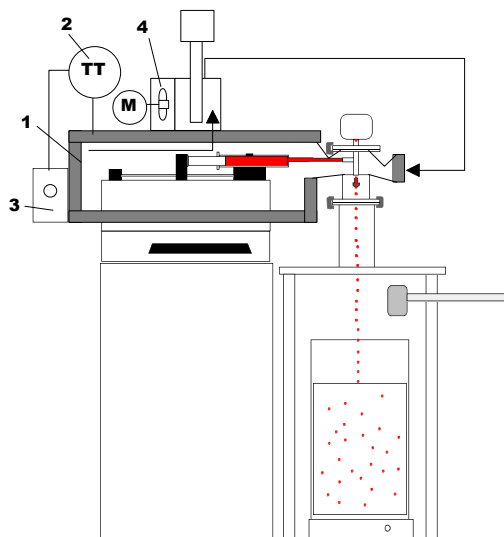

Legend:

- 1 Heating chamber (can be opened from the top to insert the syringes)
- 2 Temperature sensor inside the chamber
- 3 Temperature controller
- 4 Module with ventilator and electrical heating unit

Accuracy	<+-1	%
Recommended syringe sizes	30. 140	ml
N° of syringes	1. 2	
Volume / forward feed		variable
Linear forces	18	kg
Size l x w x h	230x280x140	mm
Weight	4.0	kg
Electrical connection	110 and 230	V